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Application Note  
Automated Microwave-Enhanced Total Synthesis of Proteins

Page 1 of 6  
ap0239v2

## Automated Microwave-Enhanced Total Synthesis of Proteins

### Summary

Proteins and long peptides can be synthesized rapidly and efficiently by microwave-enhanced SPPS on the Liberty Blue 2.0 and Liberty PRIME™ 2.0. The use of optimized microwave SPPS incorporating a new headspace flushing technology allows for higher purity synthesis of protein sequences. This technology was demonstrated on a series of biologically relevant proteins (ubiquitin, barstar, proinsulin, collagen, I protease, and MDM2) from 76-127 amino acids which were obtained in good purity through stepwise assembly without ligation steps. High purity samples were isolated from the material by elevated temperature chromatography at 60 °C on the Prodigy™ Preparative HPLC Peptide Purification system.

### Introduction

Proteins and long peptides are critical components of biological systems and comprise many important therapeutics, but research is slowed by time-intensive expression or native chemical ligation production methods. Total synthesis by SPPS provides a direct synthetic route to target specific sequences and allows for rapid generation of analogs. However, long peptides and proteins can be challenging targets for SPPS due to the iterative accumulation of impurities and tendency for aggregation to occur. Historically, SPPS was limited to small fragments for native chemical ligation and longer sequences have been of limited accessibility.<sup>1</sup> Recently, rapid flow-based methodology has seen the significant ability to assemble long sequences in extremely quick cycle times.<sup>2</sup> However, this approach is limited by the accumulation of impurities and

Application Note  
Ultrafast Peptide Synthesis at Elevated Temperature

Page 1 of 4  
ap0131v2

## Ultrafast Peptide Synthesis at Elevated Temperature

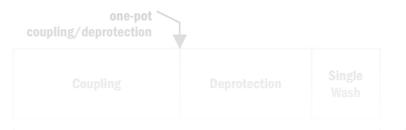
### Summary

The Liberty PRIME™ 2.0 microwave peptide synthesizer is the most advanced system available for peptide synthesis. It is based on the use of a new one-pot coupling and deprotection methodology for solid phase peptide synthesis that drastically reduces cycle time and waste usage to unprecedented levels. The system provides a complete cycle time of only 2 min 10 sec (for all 20 standard Fmoc amino acids) with only 8 mL of chemical waste produced, half the amount of CEM's highly efficient Liberty Blue 2.0 peptide synthesizer. The Liberty PRIME 2.0 HT24 is an ideal system for high-throughput peptide synthesis of both standard and complex peptides, with the ability to automatically synthesize up to 24 peptides in a day.

### Introduction

Traditional solid phase peptide synthesis involves the use of iterative and expensive deprotection and coupling steps with

side reactions at the resin bound amino functionality. The Fmoc removal then proceeds uninterrupted at elevated temperature. An optimized use of reagents results in an essentially neutral reaction mixture towards the end of deprotection step. This new procedure offers several advantages such as (a) approximately 90% reduction in solvent requirement for the deprotection step, (b) 75% reduction in solvent requirement for post-deprotection washings, (c) faster deprotection step, since the microwave ramp time is not needed, and (d) shorter cycle time due to absence of post-coupling drain step.

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